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
-

Líneas base

Versión 5.30

Multi-species

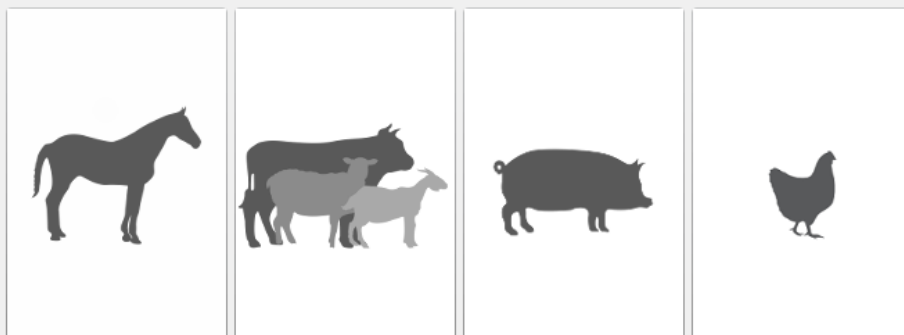ID.SOFT

With you at every step

Propósito de este manual:

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Baselines X

Calculate a baseline

Select the disease, the age of the flocks and their type (broilers, breeders, etc.) to calculate the baseline.

Label:

From: to:

Disease: Type of vaccine:

Flocks age: Between and Flock type: Elements to print: Chart List

Quick search

	Flock code	Age	Customer	Name of farm	House
	2	35 Day(s)	Client A	Client A	
	1	36 Day(s)	Client A	Client A	
	3	36 Day(s)	Client A	Client A	
	4	35 Day(s)	Client A	Client A	
	6	36 Day(s)	Client A	Client A	
	7	36 Day(s)	Client A	Client A	
	8	36 Day(s)	Client A	Client A	
	9	36 Day(s)	Client A	Client A	
	10	36 Day(s)	Client A	Client A	
	11	36 Day(s)	Client A	Client A	
	14	36 Day(s)	Client A	Client A	

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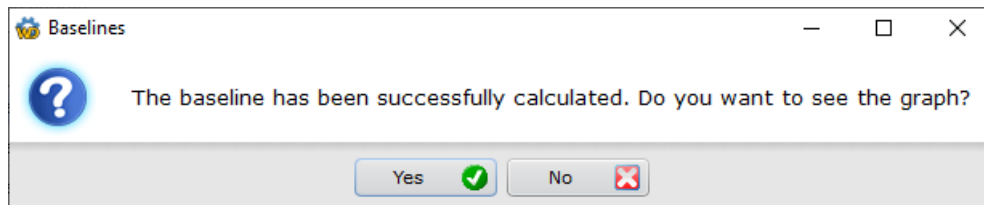
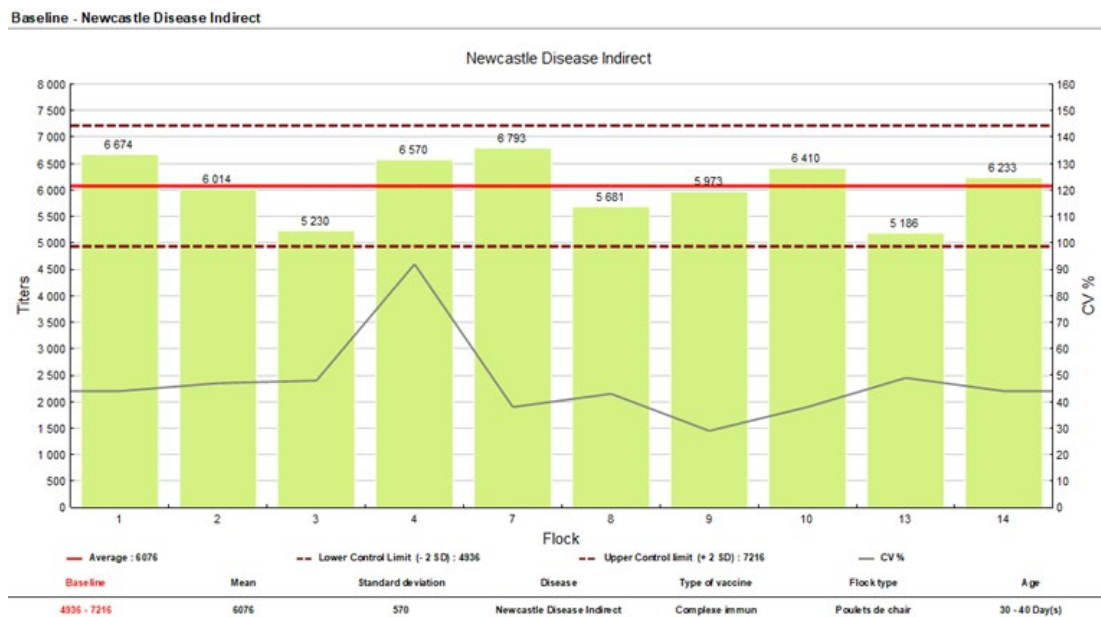
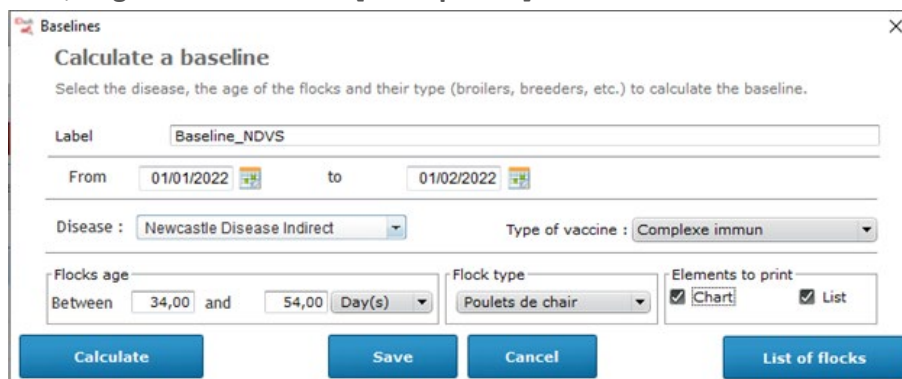
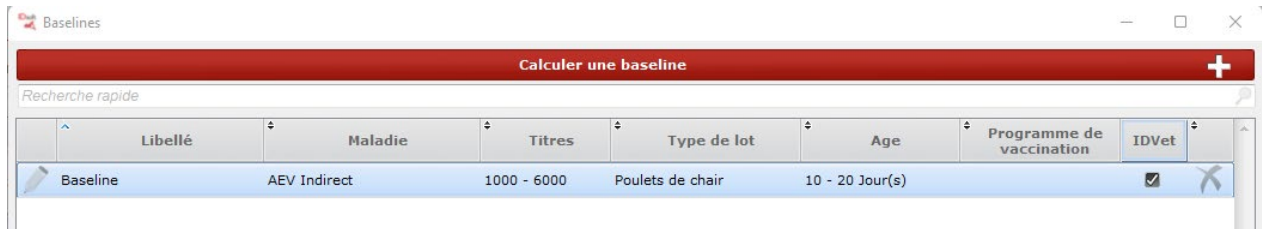
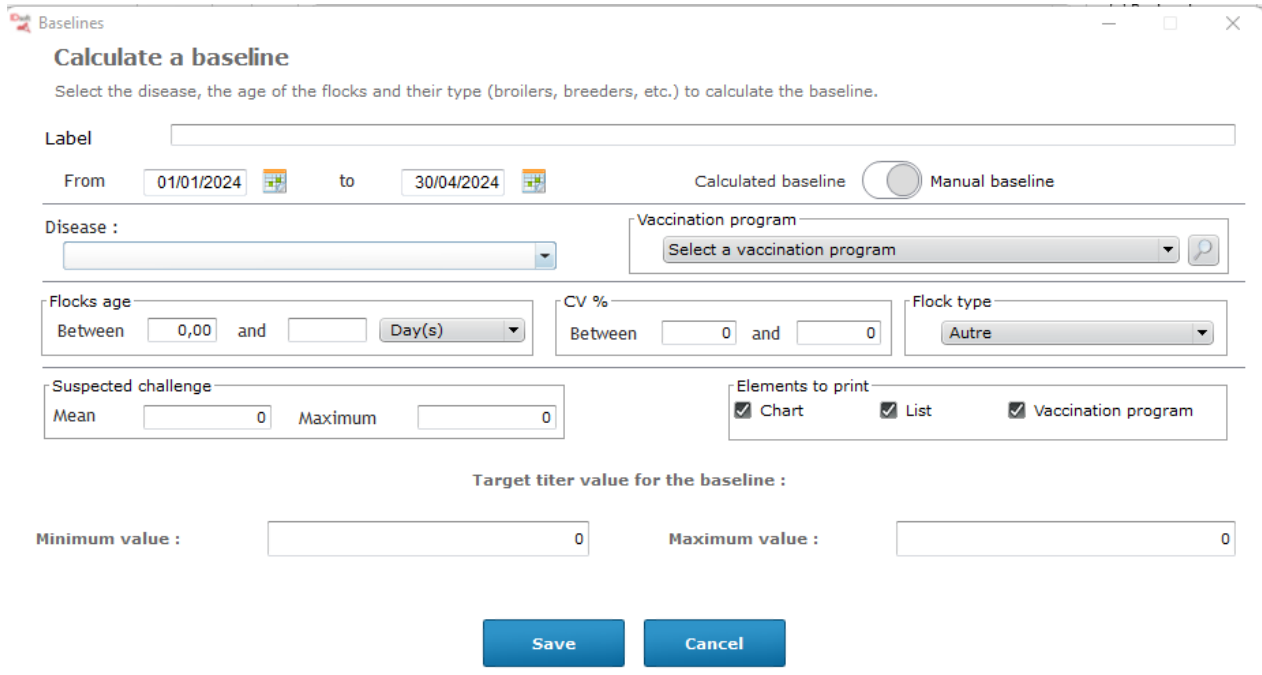
A screenshot of a form titled 'Calculate a baseline'. The form includes a 'Label' input field. Below it are 'From' and 'to' date pickers with values '01/01/2024' and '30/04/2024' respectively. There are radio buttons for 'Calculated baseline' and 'Manual baseline', with 'Manual baseline' selected. The 'Disease' field is a dropdown menu. The 'Vaccination program' field is a dropdown menu with a search icon. There are three sections for 'Flocks age', 'CV %', and 'Flock type', each with 'Between' and 'and' input fields and a unit dropdown. The 'Flocks age' unit is 'Day(s)'. The 'CV %' unit is '%'. The 'Flock type' dropdown has 'Autre' selected. There are 'Suspected challenge' fields for 'Mean' and 'Maximum', both with '0' entered. The 'Elements to print' section has checkboxes for 'Chart', 'List', and 'Vaccination program', all of which are checked. At the bottom, there are 'Minimum value' and 'Maximum value' input fields, both with '0' entered. Finally, there are 'Save' and 'Cancel' buttons.

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