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
-

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Versión 5.30

Multi-species

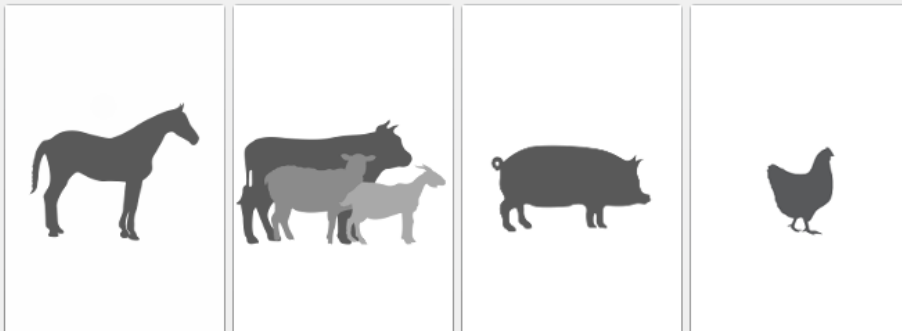ID.SOFT

With you at every step

Propósito de este manual:

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The 'Customer file' window contains the following fields and sections:

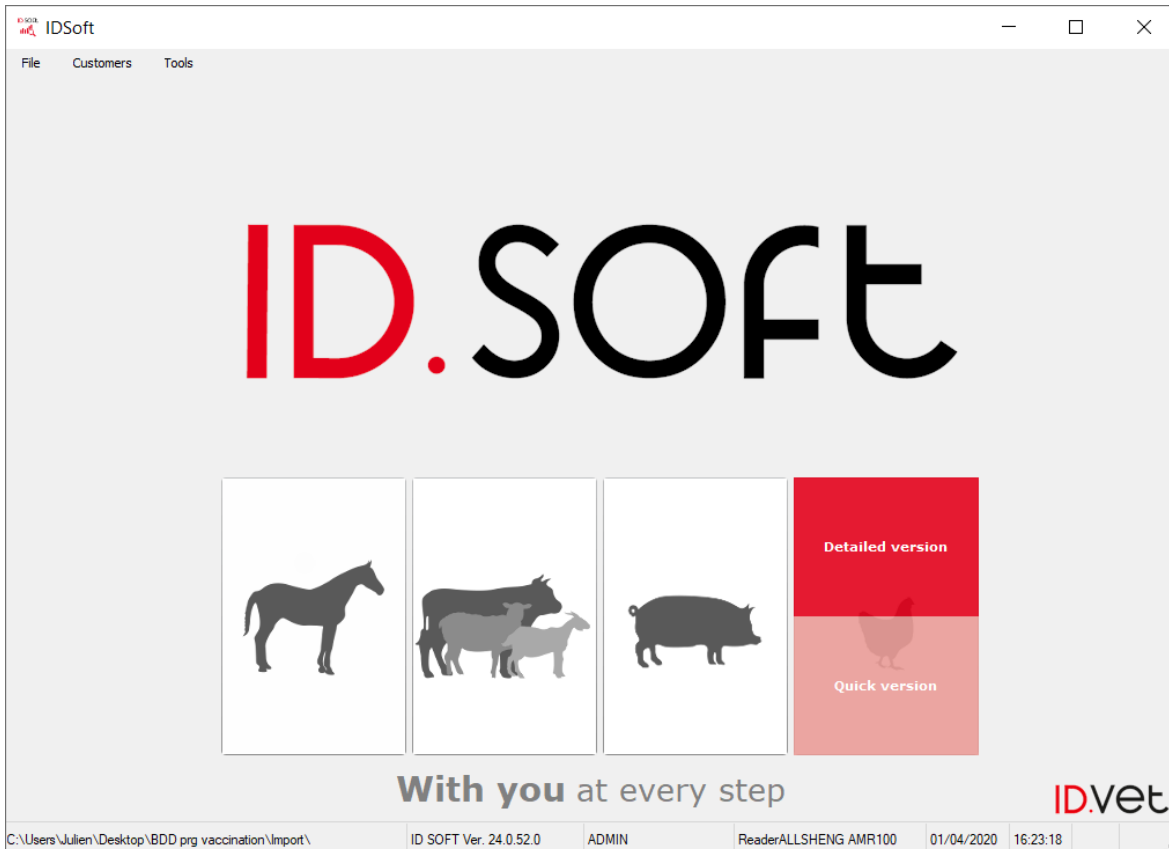
- Customer type: Breeder (dropdown), Veterinarian (checkbox), No veterinarian (checkbox)
- Name of farm: [text input]
- Name of customer: [text input]
- Address: [text input]
- Postal code: [text input], City: [text input]
- Country: [text input]
- Phone: [text input], Fax: [text input]
- Email: [text input]
- Bar code: [text input]
- Bottom section: Name of farm (selected), Flocks, Tests
- Table: Add name of farm (+)
 - Columns: Name of farm, Farm No, Location
 - Rows: 5 empty rows for data entry
- Buttons: Delete all files, Done, Cancel

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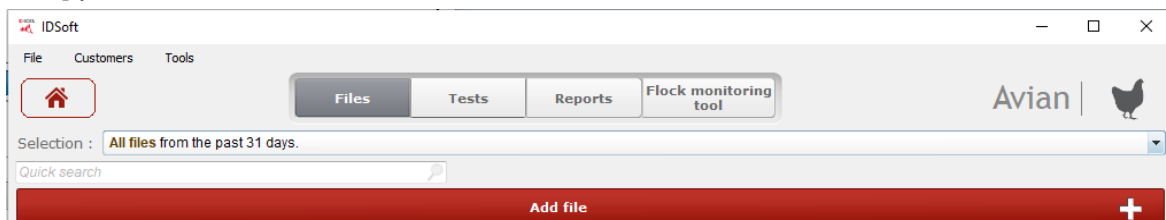
The screenshot shows a software window titled "File" with the following fields and controls:

- Date: 01/04/2020, Hour: 16:28, Operator: ADMIN
- * File No: 20200401-18
- Avian
- * Other: FARM A (with search icon), Customer site: FARM A
- Veterinarian: No veterinarian (with search icon)
- Information / Further information:
 - Name of farm: FARM A (with search icon)
 - House: 4 (with search icon)
 - Flock code: Select a flock number (with search icon)
 - Flock type: [empty]
 - Chick Hatch date: 30/03/2020
 - Bleeding date: 01/04/2020
 - Flock age: 3.00 Day(s)
- Column display: * Sample count: 0 (with green checkmark)
- Add Diseases (with plus icon)
- Table with columns: ID, Bar code, Sex
- * Required fields
- Done, Cancel

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Information Further information

Name of farm: FARM A House: 4

Flock code: Select a flock number Flock type:

Chick Hatch date: 30/03/2020 Bleeding date: 01/04/2020 Flock age: 3.00 Day(s)

Column display * Sample count: 10 Add Diseases +

ID	Bar code	Sex
01		
02		
03		
04		
05		
06		

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Column display * Sample count: 20

ID	Bar code	Sex
01	BC0001	Female
02	BC0002	Male
03	BC0003	Male
04	BC0004	Female
05	BC0005	Female
06	BC0006	Female

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ID Screen® IBD Indirect

Indirect ELISA for the detection of anti-IBD antibodies in chicken sera.

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The screenshot shows the 'Avian' software interface for file management. At the top, it displays 'File n° 20200401-18 (Pending files)' and the operator 'ADMIN'. The date is set to '01/04/2020' and the time to '16:28'. The file number '20200401-18' is highlighted in red. The interface is divided into several sections: 'Other' with a search for 'FARM A', 'Veterinarian' with 'No veterinarian', and 'Information' with fields for farm name, flock code, hatch date, house, flock type, bleeding date, and flock age. A table below shows a list of samples with IDs 01 through 06, each with a red 'X' icon. At the bottom, there are three buttons: 'Done', 'Cancel', and 'Duplicate'.

ID	Bar code	Sex
01		
02		
03		
04		
05		
06		