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
-

## Reportes

Versión 5.30

Multi-species
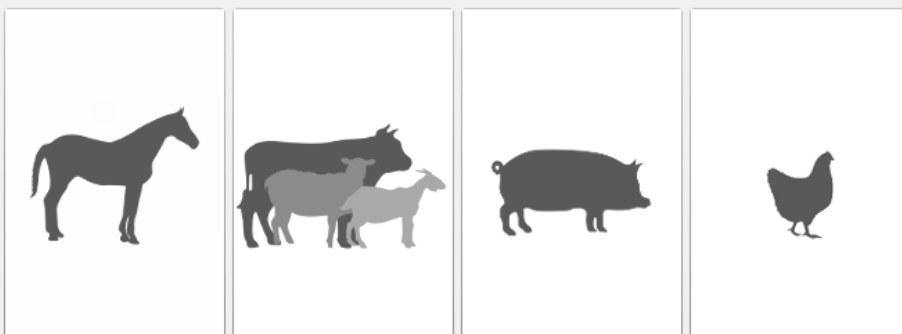
# ID.SOFT

**With you** at every step

### Propósito de este manual:

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Signature

Signature date

Species	<b>Avian</b>	Breed	<input type="text"/>
Name of farm	<b>FARM B</b>	House	<b>7</b>
Flock code	<input type="text"/>	Type	<input type="text"/>
Application	<b>Serum / Plasma - All applications</b>	Flock age	<b>3 Day(s)</b>
Test purpose	<input type="text"/>	Bleeding date	<b>08/03/2020</b>

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**EDS Indirect**

File List Graph

Serum / Plasma - All applications **Negative** <=0.5< Positive

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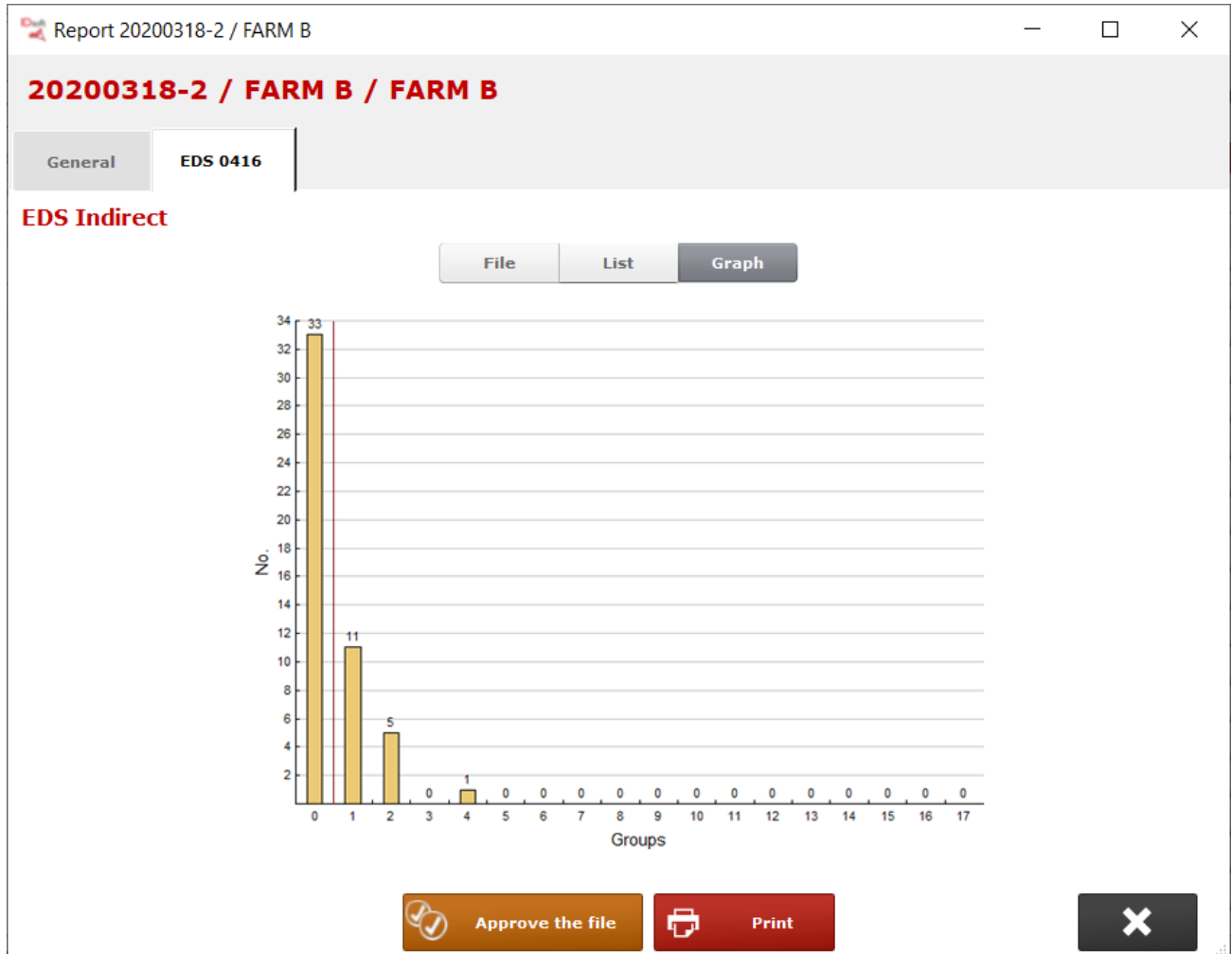
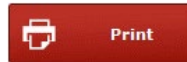
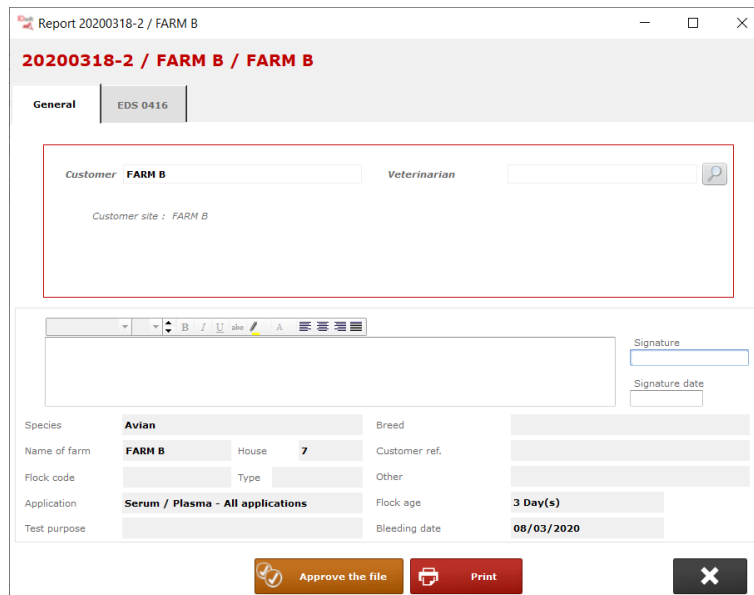
Signature date: [ ]

Species: Avian Breed: [ ]

Name of farm: FARM B House: 7 Customer ref.: [ ]

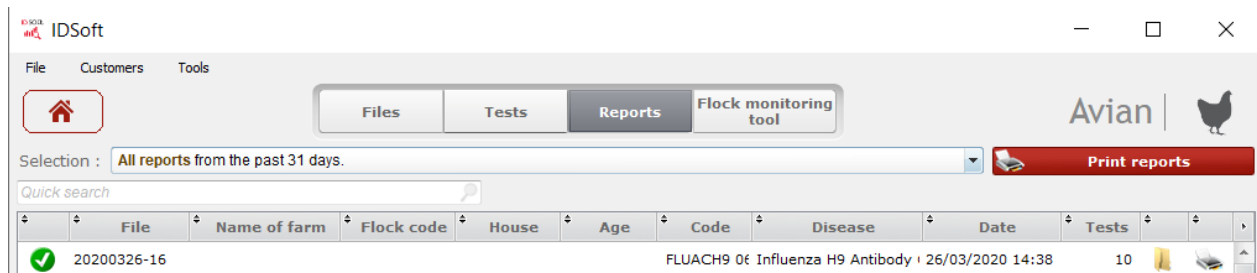
Flock code: [ ] Type: [ ] Other: [ ]

Application: Serum / Plasma - All applications Flock age: 3 Day(s)

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IDSoft Avian

File Customers Tools

Files Tests Reports Flock monitoring tool

Selection: All reports from the past 31 days. Print reports


Quick search

File	Name of farm	Flock code	House	Age	Code	Disease	Date	Tests
20200326-16						FLUACH9 0€ Influenza H9 Antibody	26/03/2020 14:38	10

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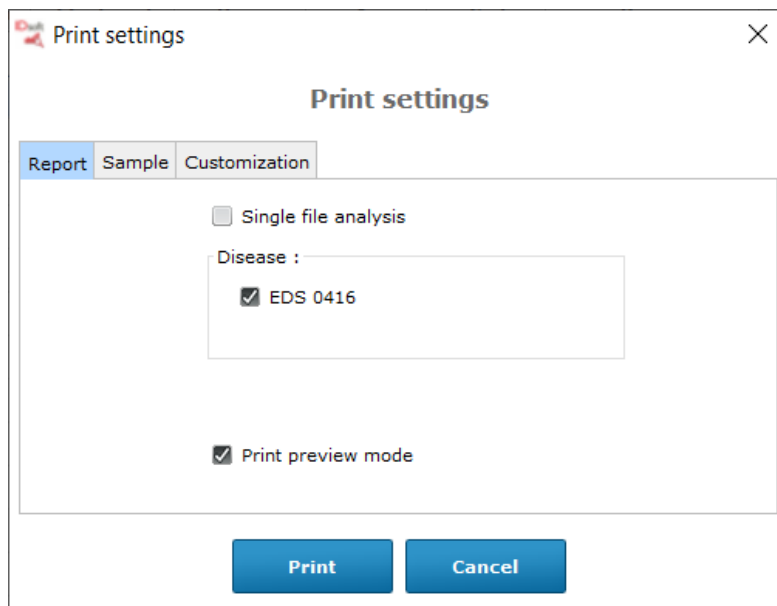
The screenshot shows a 'Print settings' dialog box with a close button (X) in the top right corner. The title bar reads 'Print settings'. Below the title bar, there are three tabs: 'Report' (selected), 'Sample', and 'Customization'. The main content area contains the following options:

- Single file analysis
- Disease :
  - EDS 0416
- Print preview mode

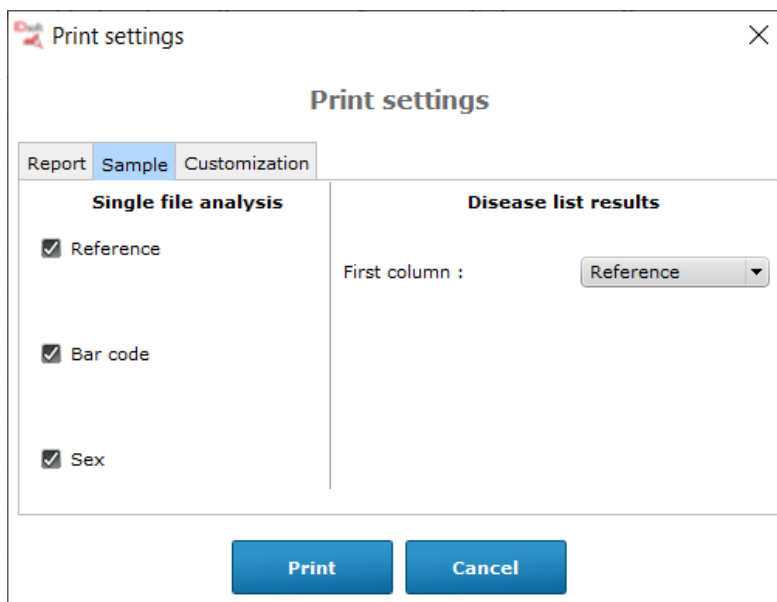
At the bottom of the dialog box, there are two blue buttons: 'Print' and 'Cancel'.

The screenshot shows the same 'Print settings' dialog box, but with the 'Sample' tab selected. The main content area is divided into two sections:

- Single file analysis**
  - Reference
  - Bar code
  - Sex
- Disease list results**
  - First column :

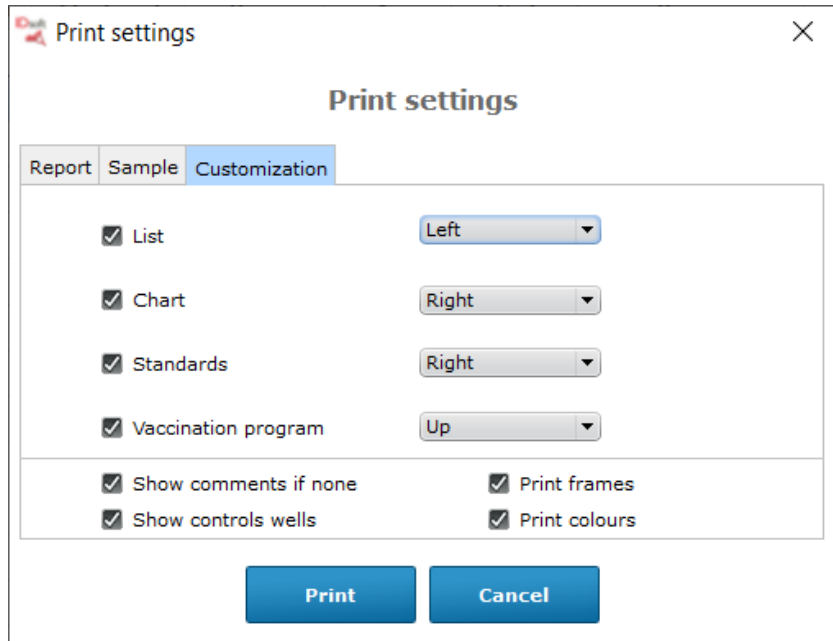At the bottom of the dialog box, there are two blue buttons: 'Print' and 'Cancel'.

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